

CHEMTEQ[®]

DG Water Test

Hydrazine

Ultra-Low and Low Range Wide Span
(PN: 83301)



Manual

CHEMTEQ[®]

600 West 24th Street, Suite B
Norfolk, Virginia 23517, USA
Tel: 757-622-2223
Toll-free: 855-CHEMTEQ (855-243-6837)
sales@chemteq.net
www.chemteq.net



1. Application

The DG Water Test PN: 83301 is a colorimetric water test kit designed to provide real-time indication and quantification of hydrazine in water.

2. Components of the Test Kit

- a. 30 Sample Vials, PN: 83301-1030
- b. One Developing Solution Bottle, PN: 83301-2000
- c. 30 Disposable Pipettes, PN: 83301-3030
- d. One Color Comparator, PN: 83301-6000

3. Components of the Test Kit Refill

- a. 30 Sample Vials, PN: 83301-1030
 - b. One Developing Solution Bottle, PN: 83301-2000
 - c. 30 Disposable Pipettes, PN: 83301-3030
- To Reorder: PN: 83301-5000

4. Specifications

4.1. Overall Specification

A. Sample Vial

- a. Weight: 0.75g (0.027oz)
- b. Dimensions: 3.3cm (1.3in), Φ : 10mm (0.4" in)
- c. Operating temperature: 4°C to 60°C (39°F to 140°F)
- d. Minimum detectable limit: 0.002ppm in 5 minutes
- e. Detection range with color comparator: 0.002 to 500ppm
- f. Color change: Colorless to yellow or orange
- g. Storage temperature: 4°C to 25°C, (39°F to 77°F)
- h. Shelf life: 1 year

B. Developing Solution

- a. Weight: 9.g (0.32oz)
- b. Dimensions: 2.3cm (0.9in), Φ : 18mm (0.69" in)

C. Color Comparator

- a. Weight: 28g (1oz)
- b. Dimensions: 10.9cm (4.3in) x 10.5cm (4.1in) x 0.7cm (0.3in)
- c. Range: 0.002ppm to 500ppm
- d. Incruments (ppm): 0.002, 0.005, 0.01, 0.02, 0.03, 0.04, 0.05, 0.06, 0.07, 0.08, 0.09, 0.1, 0.2, 0.3, 0.4, 0.5, 0.75, 1.0, 2.0, 4.0, 6.0, 8.0, 10.0, 15, 25, 50, 75, 100, 125, 150, 175, 200, 225, 250, 275, 300, 325, 350, 375, 400, 425, 450, 500.

D. Disposable Pipette

- a. Weight: 0.75g (0.027oz)
- b. Dimensions: 3.4cm (2.5in), Φ : 6.8mm (0.27" in)
- c. Capacity: 0.2ml

4.2. Cross interferences

Aromatic amines and aromatic hydrazine produce similar color. No other interferences known

5. Operating Instructions

- a. Use gloves and protective glasses when handling Hydrazine.
- b. Ensure that packaging pouch is intact.
- c. Open packaging pouch by tearing off the top part from one of side notches.
- d. Remove one sample vial from packaging pouch.
- e. Open sample vial cap and add 0.4ml of water sample using the disposable pipette (Figure 1).
- f. Add three drops of developing solution to sample vial (Figure 2).
- g. Firmly close sample vial cap and shake (Figure 3).
- h. Wait 5 minutes with occasional shaking for complete color development.

NOTE: Reading the result before 5 minutes or waiting much longer than 5 minutes may lead to negative or positive results

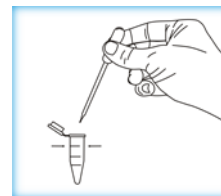


Figure 1



Figure 2

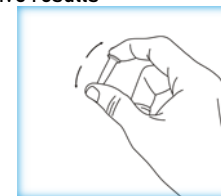


Figure 3



Figure 4

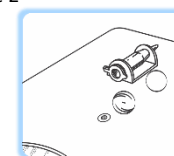


Figure 5



Figure 6

- i. Use the color comparator (Figure 6) to determine the concentration of hexavalent chromium in the water sample
 - a. Insert the sample vial into the vial holder as shown in Figures 4 and 5.
 - b. Turn the bottom color wheel to match colors. The color formed in the sample vial is directly proportional to the concentration of hexavalent chromium in the water sample.
 - c. If the color of the sample vial is darker than the darkest color on the bottom wheel. Reset the bottom wheel to the black diamond and turn top wheel anticlockwise until color match obtained.
 - d. If the color of the sample vial still is darker than the darkest color of the top wheel, dilute your sample with equal amount of DI water and repeat steps (a) to (c), make sure to multiply the reading on the wheel by two after diluting your sample