


DG Filter 700 Series with BTI 2 Use and Refilling Instructions



DG Filter 700 S Use & Refilling Instructions

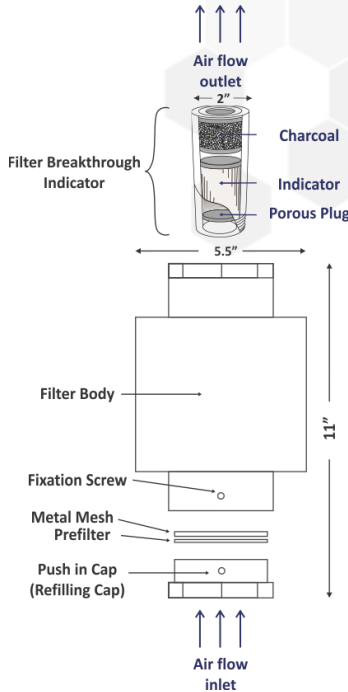

Instruction of Use

1. Remove the DG Filter from packaging bag.
2. Wipe the unit with a damp cloth.
3. Plug into a grounded 110-volt outlet, run on TURBO (high) for a few minutes and then choose the desired setting.
4. Always let the DG Filter run for an hour or more before using in the desired area.
5. Connect suitable Breakthrough Indicator (BTI) to the filter outlet, refer to BTI manual for details
6. Replace or refill filter when BTI change color.

Filter Refilling Instructions

1. Remove fixation screw.
2. Remove refilling cap, prefilter and sorbent retainer (metal mesh).
3. Empty consumed sorbent disposes of in accordance with all applicable regulations.
4. To charge the empty filter with fresh sorbent media:
 - a. Choose the suitable sorbent media
 - b. Charge the filter with about one pound of fresh sorbent and shake the filter with occasional pressing to ensure uniform settling of sorbent in the filter.
 - c. Repeat step (b) three times.
 - d. Replace sorbent retainer and prefilter.
 - e. Align the fixation screw opening on the refilling cap with the opening on the filter.
 - f. Push in the refilling cap and screw in the fixation screw.

CHEMTEQ[®]

Chemteq, Inc., 600 West 24th Street, Suite B, Norfolk, Virginia 23517, Tel 757-622-2223, sales@chemteq.net
©2009 Chemteq, Inc.

MP-7219000-092420-C